

FireSting[®]-PRO

Optical Multi-Analyte Meter

USER MANUAL

O₂

pH

T



FireSting®-PRO

Optical Multi-Analyte Meter

Document Version 1.12

The FireSting®-PRO is released by:

PyroScience GmbH

Kackertstrasse 11
52072 Aachen
Germany

Phone +49 (0)241 5183 2210
Fax +49 (0)241 5183 2299
Email info@pyroscience.com
Web www.pyroscience.com

Registered: Aachen HRB 17329, Germany


TABLE OF CONTENT

1	Overview	4
2	Introduction	5
3	Quick Start	6
4	Connecting Sensors	7
5	Optical Sensors	8
	5.1 Fiber-optic Sensors.....	8
	5.2 Contactless Sensors	9
6	Extension Port	11
	6.1 Connector X1 (Power, Digital Interface, Analog In)	11
	6.2 Connector X2 (Analog Output)	12
7	Specifications	14
	7.1 Mounting Hole Dimensions of the FireSting® meters	15
8	Related Documents	16
9	Warnings & Safety Guidelines	17

1 OVERVIEW

The compact USB-powered fiber-optic meter **FireSting®-PRO** with 1, 2, or 4 channels for multiple analytes and sensor heads features:

- freely configurable channels for O₂, pH and temperature
 - broad optical sensor portfolio (multiple fiber-based and contactless sensor heads for several ranges)
 - with (ultra-)high speed sampling
 - zero-noise and zero-latency temperature compensation
 - improved ambient light suppression and
 - smart measuring modes for prolonged sensor lifetime
 - pH, oxygen and temperature determination simultaneously in one sample
-

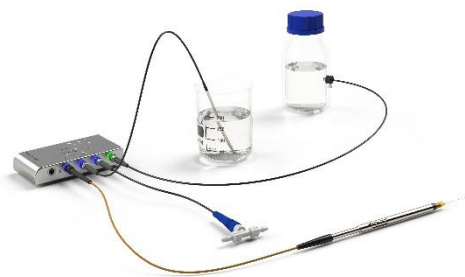

This single-device solution for multi-analyte measurements has integrated atmospheric pressure and humidity sensors for precise and easy oxygen sensor calibration, but also for automatic pressure compensation of the oxygen measurements. Furthermore, the **FireSting®-PRO** offers 4 analog outputs and a broadcast-mode.


The new innovative and user-friendly **Pyro Workbench** allows operation of several **FireSting®-PRO** meters in parallel as a scalable multi-channel system.

2 INTRODUCTION

The FireSting®-PRO comes with 1, 2, or 4 channels (optical sensor ST-connectors **1** to **4**) for up to 4 fiber-optic sensors and one connector (**T**) for an external Pt100 temperature probe. The optical channel connectors are color coded, indicating the analyte (**oxygen**, (pre-calibrated) **pH** or **temperature**), which is currently measured and can be changed for each channel. The air inlet equilibrates the internal temperature, pressure and humidity sensors with the surrounding. Avoid covering these holes to ensure free air circulation towards the internal sensors.



The micro-USB connector on the left side panel provides energy supply and data exchange with the PC. Right-hand side of it, a connector **X1** for power and digital interface (7-pins) and a connector **X2** for analog output (5 pins) is located.

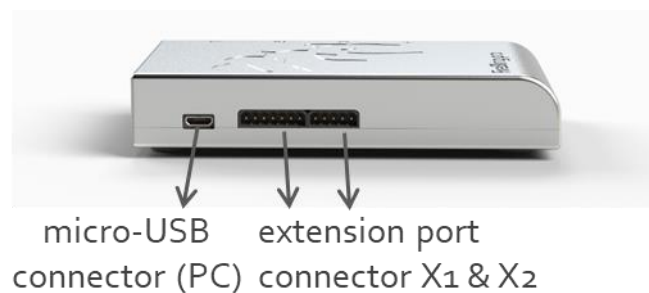

3 QUICK START

Step 1: Download the correct software and manual from the downloads tab of your purchased device on www.pyroscience.com

Step 2: Connect the **FireSting®-PRO** meter with the micro-USB cable to the Windows PC/laptop (Windows 7, 8, 10).

Step 3: Connect appropriate **PyroScience** sensor(s) to the optical sensor connectors **1** to **4** at the device (see 4).

Step 4: Connect an external temperature sensor (item no. **TDIP15** or **TSUB21**) to the Pt100 connector or, alternatively, an optical temperature sensor (see 5) to one of the optical sensor connectors **1** to **4** for automatic temperature compensation.

Step 5: Prepare appropriate calibration standards, as described in the respective sensor manuals (see 8).

Step 6: Start the logger software by clicking on the short-cut "**Pyro Workbench**" on your desktop.

Step 7: Open the **Settings Wizard** by clicking on the **FireSting®-PRO** picture. Select the **respective analyte** and enter all sensor settings for each sensor, including an appropriate mode of temperature compensation.

Step 8: Open the **Calibration Wizard** and follow the calibration instructions for each sensor. Measurements with the respective sensors will start automatically after all required sensor calibrations have been performed.

Step 9: Configure the **Graphs** according to your preferences.

Step 10: Activate **Data Logging**.

4 CONNECTING SENSORS

The fiber-optic oxygen, pH and temperature sensors, as well as optical fibers needed for read-out of contactless sensors are connected to the ST-connectors of the **FireSting®-PRO (1 to 4)** with a male fiber plug.

- Remove the black caps from the plug of the sensor / fiber.
- Remove the red caps from the sensor ports at the **FireSting®-PRO** (the red caps should be put on again if it is not in use anymore to protect the optics).
- Insert the male fiber plug of the sensor cable into the ST-port (female fiber connector) of the **FireSting®-PRO** and turn the bayonet coupling gently clockwise until the plug is locked firmly.



5 OPTICAL SENSORS

The FireSting®-PRO is compatible with a broad range of optical oxygen, pH and temperature sensors from **PyroScience**. For an overview of available optical sensor types, please see the **PyroScience** website. FireSting®-PRO devices labeled with “ready for precal pH” and all devices with SN>23360000 can be used for calibration free measurements with precalibrated pH sensors.



5.1 Fiber-optic Sensors


Sensor	Item	Analyte	Application
Miniprobes	OXROB...	O ₂	stirred water, gas
	⁺ PHROB ...	pH	water & media
Retractable Tip Minisensors	*OXR...	O ₂	water, gas & semi-solid samples
	TPR...	Temp.	
Fixed Tip Minisensors	*OXF...	O ₂	water, gas & semi-solid samples (esp. seawater)
	TPF...	Temp.	
	OXF...-PT	O ₂	gas (puncturing septa/packaging)
	PHF-PK...	pH	water & media
Bare Fiber Minisensors	OXB...	O ₂	water, gas & custom
	TPB...	Temp.	
Solvent-Resistant Probes	OXSOLV	O ₂	approved polar and non-polar solvents
	OXSOLV-PTS	O ₂	approved solvent vapors

*also as Microsensor; ⁺ different pH ranges; **water**=water, seawater, aqueous solutions


5.2 Contactless Sensors

Sensor	Item	Analyte	Application
Nanoprobes	OXNANO	O ₂	aqueous solutions & microfluidics
Sensor Spots	OXSP5	O ₂	water & gas
	TPSP5	Temp.	
	*PHSP5...	pH	water & media
Sensor Vials	OXVIAL...	O ₂	water & gas
	TOVIAL...	Temp. & O ₂	
	+PHVIAL...	pH	water & media
	+PHTVIAL...	Temp. & pH	
	+PHTOVIAL...	Temp. & pH & O ₂	
Flow-Through Cells	OXFLOW...	O ₂	water & gas
	OXFTC...		
	TPFLOW	Temp.	
	TOFTC2	Temp. & O ₂	
	+PHFLOW...	pH	water & media

⁺ different pH ranges; **water**=water, seawater, aqueous solutions

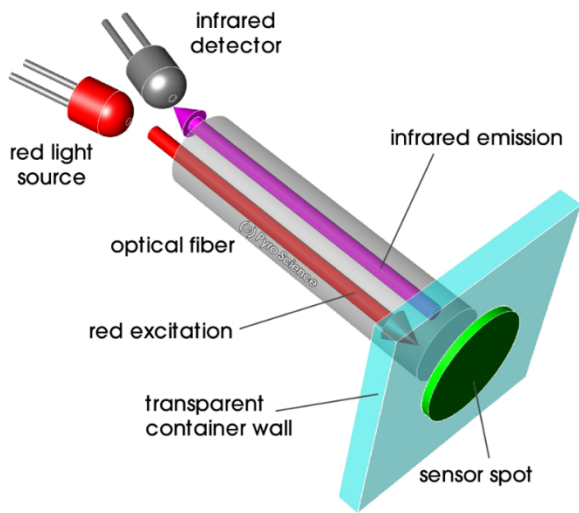

Principle of contactless read-out of sensor spots and sensor vials

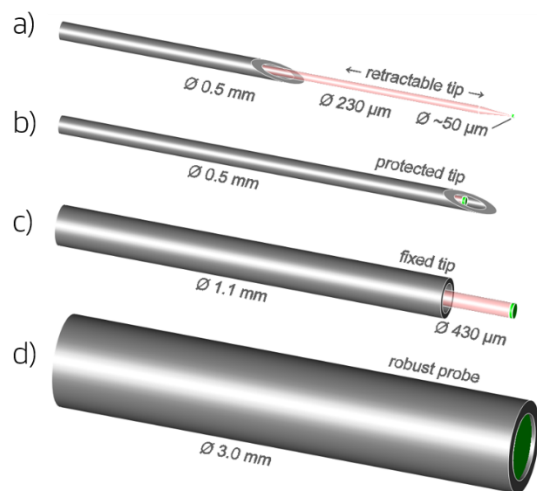


Examples of available sensor tips:

a) retractable microsensor **b)** fixed minisensor with protected tip and with **c)** protruding sensor tip, **d)** robust probe

6 EXTENSION PORT

The extension port of the **FireSting®-PRO** consists of the two connectors **X1** and **X2** (fitting connector plugs can be obtained from **Phoenix Contact** item no. **1778887** and **1778861**).

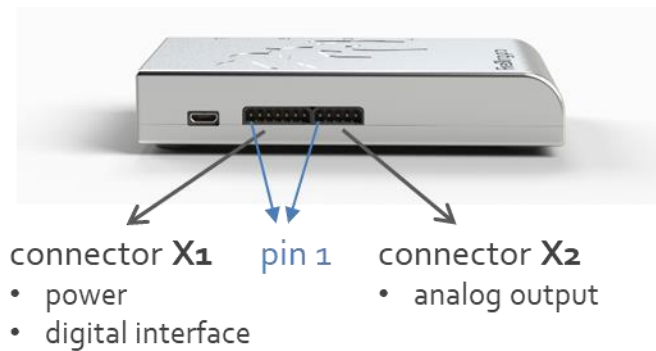

6.1 Connector X1 (Power, Digital Interface, Analog In)

The pin configuration of the connector **X1** is given in the table below. Pins 1-2 (GND and VCC) can be used for providing an external power supply (3.5...5.0 VDC), if the **FireSting®-PRO** should not be powered via the USB port. The transmit and receive pins of the UART-interface are given at pins 4 (TXD) and 5 (RXD) (communication protocol on request). When using the UART-interface it is recommended to connect pin 3 (/USB_DISABLE) to pin 1 (GND), which disables the USB interface.

Pin 6 (/PAUSE_BROADCAST) and Pin 7 (/TRIGIN) are related to the so called “broadcast mode” which can be configured in the PC control software (e.g., **Pyro Workbench** or **Pyro Developer Tool**). During broadcast mode, the device triggers itself periodic measurements which can be read out from the analog outputs or from a text message transmitted via the USB/UART interface. For more details refer to the manual of the respective control software are to the communication protocol (available on request).

Pin 7 (/TRIGIN) acts as a trigger input for the broadcast mode. Note, the option “Enable Trigin” must be enabled in the broadcast settings in the control software. Every time this pin is connected to pin 1 (GND) then an additional broadcast measurement is triggered.

Pin 6 (/PAUSE_BROADCAST) acts as a master switch of the broadcast mode. As long this pin is tied to pin 1 (GND), then the broadcast mode is paused. Neither periodic broadcast measurements, nor triggered broadcast measurements are performed.

Pin	Name	Function	Description
1	GND	Power	Ground
2	VCC	Power	Power supply, 3.5V to 5.0V DC max. 70 mA (typ. 40 mA)
3	/USB_DISABLE	Digital Input	Connect to GND for disabling the USB interface
4	TXD	Digital Output (UART TX)	UART interface with 3.3V levels (5V tolerant), baud rate 115200, 8 data bits, 1 stop bit, no parity, no handshake
5	RXD	Digital Input (UART RX)	UART interface with 3.3V levels (5V tolerant), baud rate 115200, 8 data bits, 1 stop bit, no parity, no handshake
6	/PAUSE_ BROADCAST	Digital Input (0V or 3.3V, internally pulled-up to 3.3V)	Connect to GND for pausing any broadcast-mode operation.
7	/TRIGIN	Digital Input (0V or 3.3V, internally pulled-up to 3.3V)	Triggers a broadcast measurement every time, when this pin is tied to GND.

6.2 Connector X2 (Analog Output)

The connector **X2** provides 4 analog outputs with a range of 0-2.5V DC at a resolution of 14 bits (see table below). Refer to the **Pyro Workbench** manual how to configure the analog outputs.


Pin	Name	Function	Description
1	GND		Ground
2	AO_A	Analog Output (0 - 2.5 V DC) (14-bit resolution)	Analog Output Port A (alternatively digital alarm output)
3	AO_B	Analog Output (0 - 2.5 V DC) (14-bit resolution)	Analog Output Port B (alternatively digital alarm output)
4	AO_C	Analog Output (0 - 2.5 V DC) (14-bit resolution)	Analog Output Port C (alternatively digital alarm output)
5	AO_D	Analog Output (0 - 2.5 V DC) (14-bit resolution)	Analog Output Port D (alternatively digital alarm output)

7 SPECIFICATIONS

Feature	Specification
Dimensions	68 x 120 x 22 mm (housing) 78 x 120 x 24 (total)
Weight	ca. 290 g
Interface	USB 2.0
Power supply	USB-powered (max 50mA at 5V)
Supported operating systems	Windows 7, 8, 10
Operating temperature	0 to 50°C
Max. relative humidity	non-condensing conditions
Optical sensor port	1, 2, or 4 (dependent on model)
Optical sensors	complete PyroScience sensor portfolio for O ₂ , pH & T
Optical sensor connector	fiber-optic ST-plug
Max. sample rate	ca. 10 samples per second (depending on Settings)
External temperature port Range, Resolution, Accuracy*	1 channel for 4-wire Pt100 -30°C to 150°C, 0.02°C, ±0.5°C
Internal temperature sensor Range, Resolution, Accuracy*	-40 to 125°C, 0.01°C, ±0.3°C
Internal pressure sensor Range, Resolution, Accuracy	300 to 1100 mbar, 0.1 mbar, typ. ±3 mbar
Internal humidity sensor Range, Resolution, Accuracy	0 to 100% rel. humidity (RH), 0.04% RH, typ. ±0.2% RH
Digital interface at extension port X1 (7 pins)	UART with 3.3V levels (5V tolerant), 115 200 baud, 8 data bit, 1 stop bit, no parity, no handshake

Connector plug port X1	Phoenix Contact, item no. 1778887
Analog output (4 channels) at extension port X2 (5 pins)	0 to 2.5 VDC 14-bit resolution
Connector plug port X2	Phoenix Contact, item no. 1778861

7.1 Mounting Hole Dimensions of the FireSting® meters

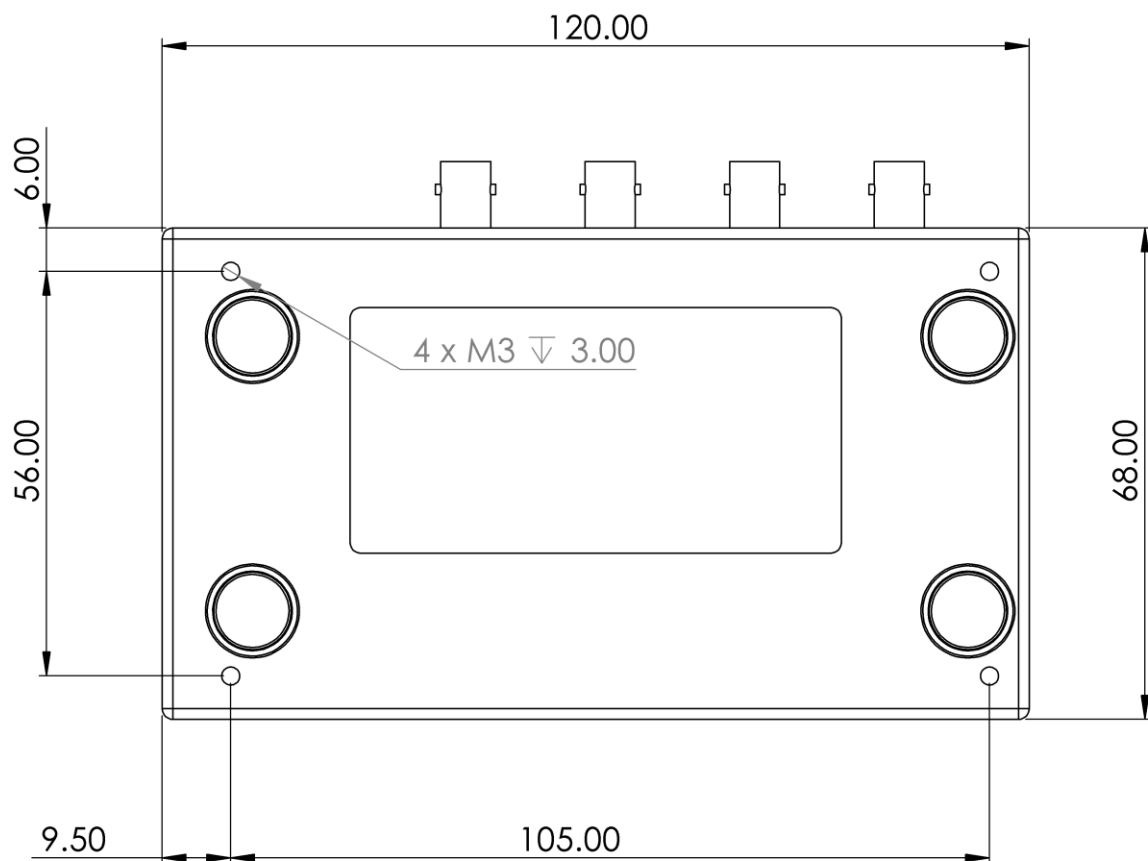

8 RELATED DOCUMENTS

Detailed instructions for using the **Pyro Workbench** and application of optical oxygen, pH and temperature sensors:

- manual for logger software "**Pyro Workbench**" (Windows)
- manuals for optical sensors from **PyroScience** (oxygen, pH, temperature)

9 WARNINGS & SAFETY GUIDELINES

Before using the FireSting®-PRO and its sensors, read carefully the instructions and user manuals.

In case of problems or damage, disconnect the instrument and mark it to prevent any further use! Consult **PyroScience** for advice! There are no serviceable parts inside the device. Please note that opening the housing will void the warranty!

The FireSting®-PRO is not watertight, is sensitive to corrosive conditions and to changes in temperature causing condensation. Avoid any condition (e.g. direct sun light) causing a heating of the device above 50°C (122°F) or below 0°C (32°F). Avoid any elevated humidity causing condensing conditions.

Handle the sensors with care especially after removal of the protective cap! Prevent mechanical stress to the fragile sensing tip! Prevent injuries with needle-type sensors!

Calibration and application of the sensors is on the user's authority, as well as data acquisition, treatment & publication!

The optical sensors and meter **FireSting®-PRO** are **not** intended for medical, diagnostic, therapeutic, or military purposes or any other safety-critical applications. The sensors must **not** be used for applications in humans and must not be brought in direct contact with foods intended for consumption by humans.

The FireSting®-PRO and optical sensors should be used in the laboratory by qualified personnel only, following the user instructions and the safety guidelines of the manual, as well as the appropriate laws and guidelines for safety in the laboratory!

Keep the sensors and the fiber-optic meter **FireSting®-PRO** out of reach of children!

CONTACT

PyroScience GmbH

Kackertstraße 11
52072 Aachen
Deutschland

Tel.: +49 (0)241 5183 2210

Fax: +49 (0)241 5183 2299

info@pyroscience.com

www.pyroscience.com